VWR

Stereomicroscope  **VisiScope®** series 300

INSTRUCTION MANUAL

Model	European Catalogue Number
SZB350	630-1578
SZT350	630-1580
SZB350 OH	630-1583
SZT350 OH	630-1584

Version: 3
Issued: 19, 05, 2014

Legal Address of Manufacturer

Europe

VWR International bvba
Researchpark Haasrode 2020
Geldenaaksebaan 464
B-3001 Leuven
+ 32 16 385011
<http://be.vwr.com>

Country of origin: **ITALY**

Table of Contents

Warning

Safety information

Package Contents

Unpacking

Intended Use

Symbols and conventions

Specifications

Instructions for use

Overview

Description of Buttons and Switches

Operation

Troubleshooting

Repair and maintenance

Accessories and spares

Technical Service

Warranty

Compliance with local laws and regulations

Disposal

Warning

This microscope is a scientific precision instrument designed to last for many years with a minimum of maintenance. It is built to high optical and mechanical standards and to withstand daily use.

We remind you that this manual contains important information on safety and maintenance, and that it must therefore be made accessible to the instrument users.

We decline any responsibility deriving from incorrect instrument use that does not comply with this manual.

Safety Information

Avoiding Electrical Shock

Before plugging in the power supply, make sure that the supplying voltage of your region matches with the operation voltage of the equipment and that the lamp switch is in off-position.

Users should observe all safety regulations of the region. The equipment has acquired the CE safety label. However, users do have full responsibility to use this equipment safely.

Please follow the guidelines below, and read this manual in its entirety to ensure safe operation of the unit.

Package Contents

DESCRIPTION	QUANTITY
Microscope stand with focus and illuminators (NO illumination on -OH versions)	1
Optical stereo head	1
Eyepiece WF10x/20mm	2
Disc for transmitted illumination (NOT included in -OH versions)	1
Dust cover	1

Unpacking


The microscope is housed in a moulded Styrofoam container. Remove the tape from the edge of the container and lift the top half of the container. Take some care to avoid that the optical items (head and eyepieces) fall out and get damaged. Using both hands (one around the arm and one around the base), lift the microscope from the container and put it on a stable desk.

Place the observation head onto the top of the arm and tighten the lock-screw. Insert the eyepieces into the eye tubes.

Connect the provided power supply cable to the input socket on the rear of the microscope.

Overhanging stand assembling


- Screw the main long pole into the thread on the base.

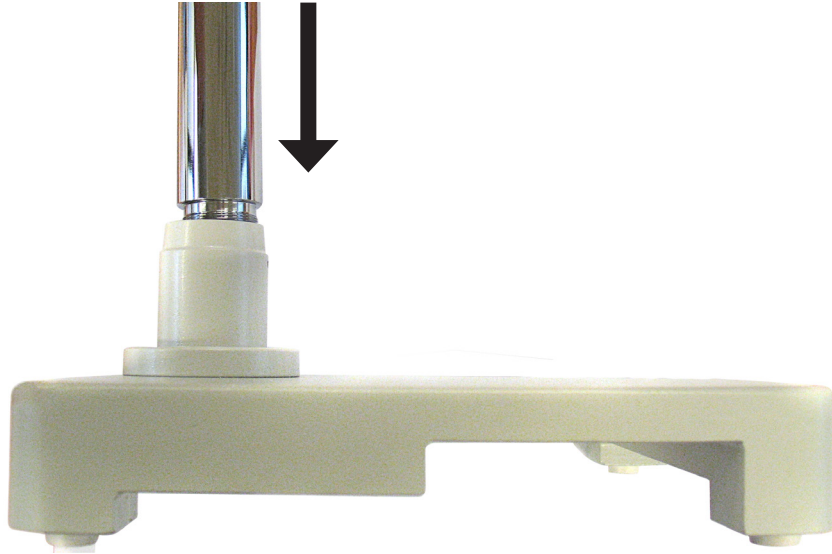
- Slide the stop-ring along the pole, and fix it with a provided screw.


- Prepare the horizontal rod, with 4 screws.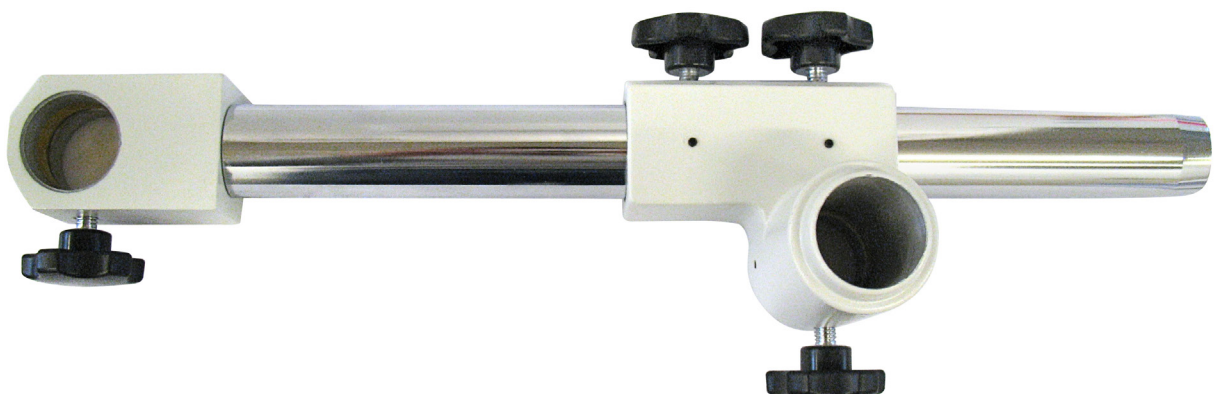


-
- Then slide the rod along the main pole. Make it rest on the stop-ring, it will provide additional safety, avoiding that the microscope could hit the sample.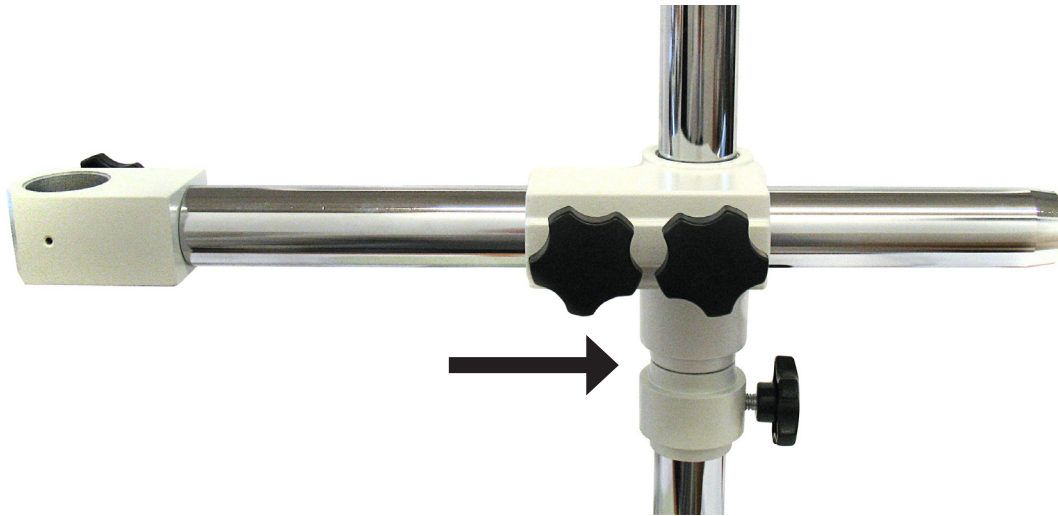


- At the end of the horizontal arm, slide the short rod in.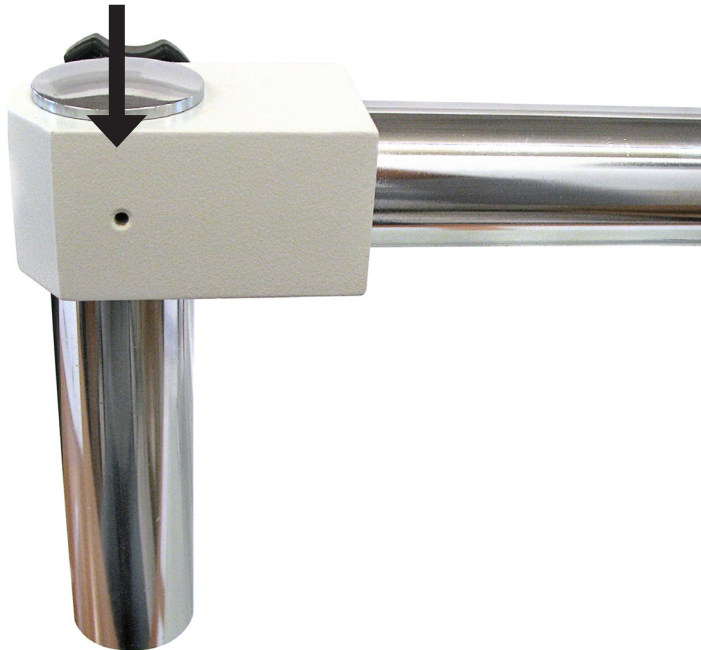

- Take the focus, and mount it on the short rod. At the bottom of the rod, screw the fixing cap.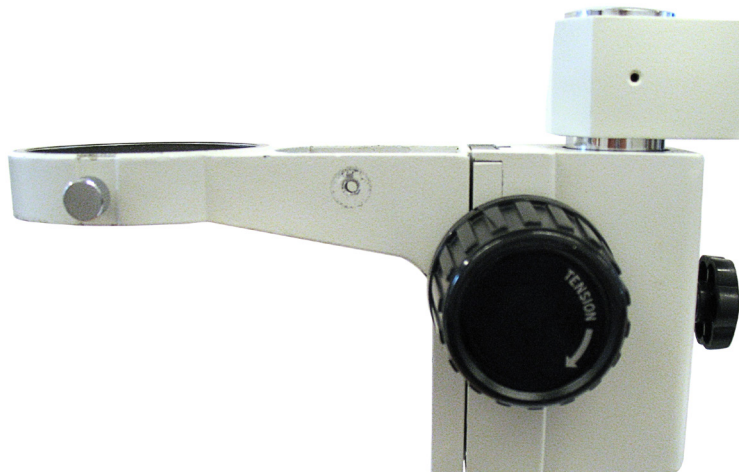

- Finally, mount the stereomicroscope head into the focus, and lock the head with the provided screw.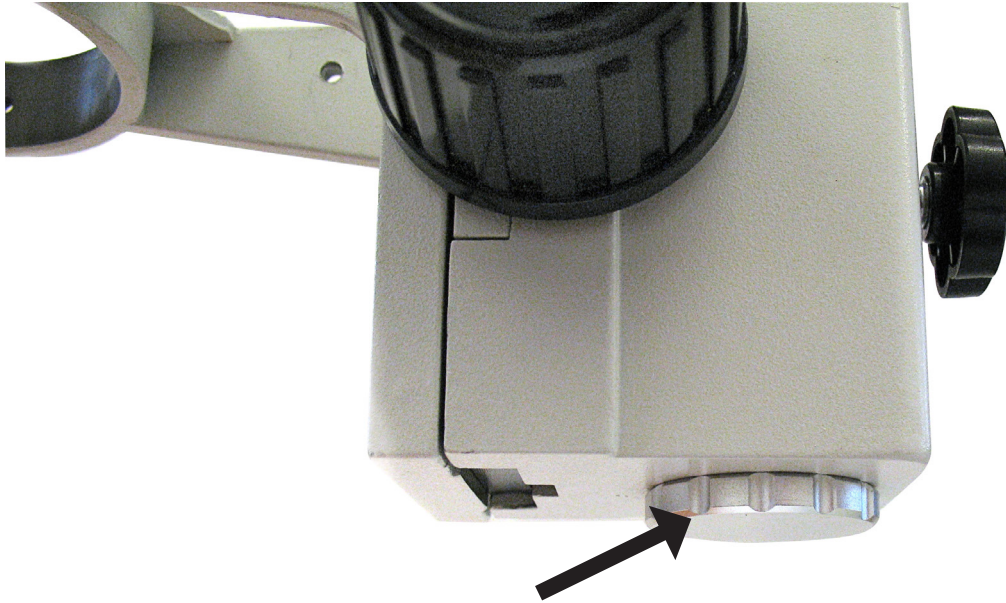

Intended use

For research use only. Not intended for any animal or human therapeutic or diagnostic use.

Symbols and conventions


The following chart is an illustrated glossary of the symbols that are used in this manual.

	CAUTION This symbol Indicates a potential risk and alerts you to proceed with caution
---	---

PRODUCT SPECIFICATIONS

Model	Head	Eyepieces	Objectives	Working Distance	Stand	Illumination
SZB350	Binocular	EWF 10x/20mm	0,7 4,5x Zoom	100 mm	Pillar stand	Incident and transmitted 12V/15W halogen
SZT350	Trinocular	EWF 10x/20mm	0,7 4,5x Zoom	100 mm	Pillar stand	Incident and transmitted 12V/15W halogen
SZB350 OH	Binocular	EWF 10x/20mm	0,7 4,5x Zoom	100 mm	Overhanging stand	---
SZT350 OH	Trinocular	EWF 10x/20mm	0,7 4,5x Zoom	100 mm	Overhanging stand	---


OVERVIEW

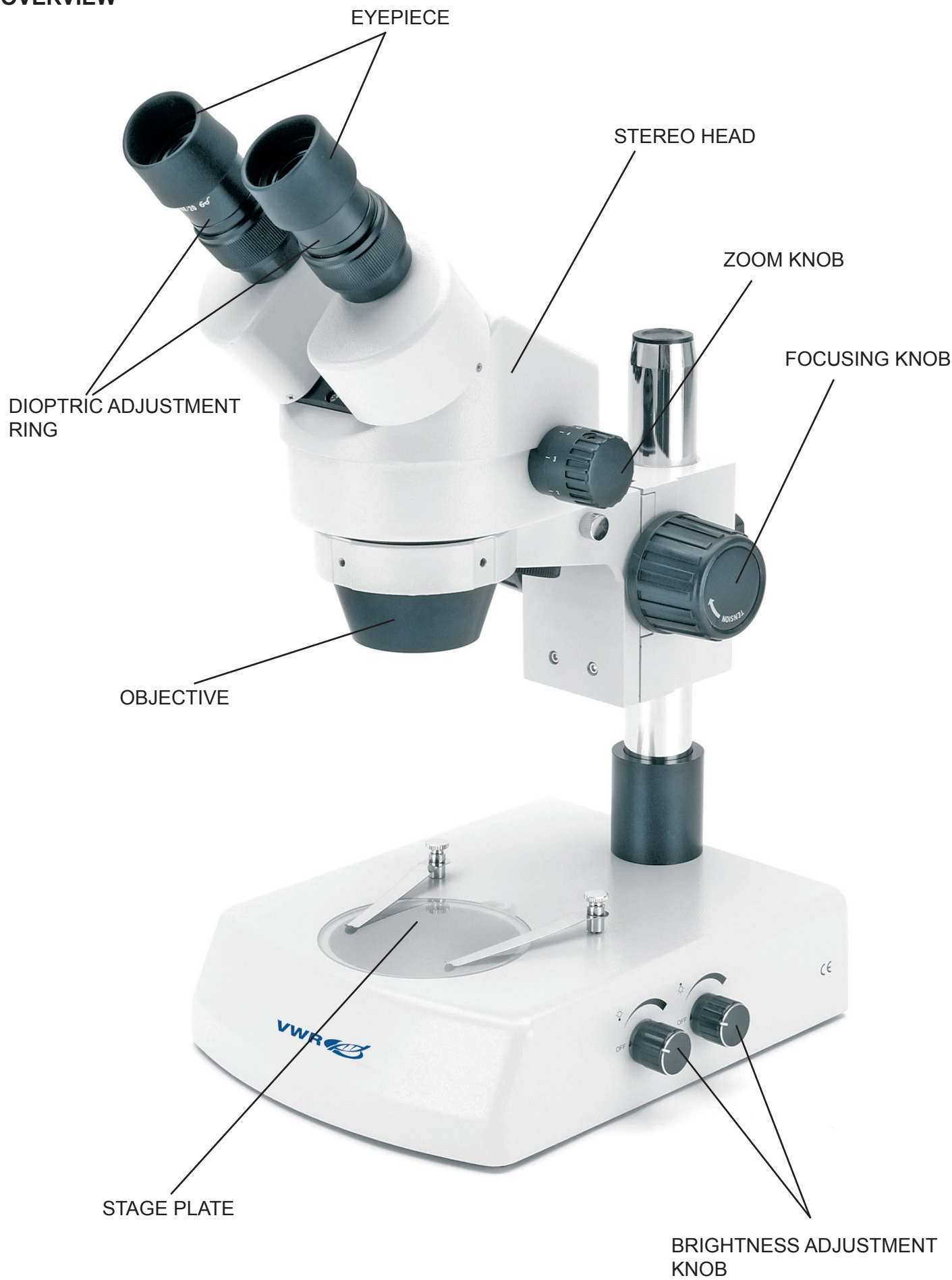Overhanging stand version (-OH)

DESCRIPTION OF BUTTONS AND SWITCHES Front View

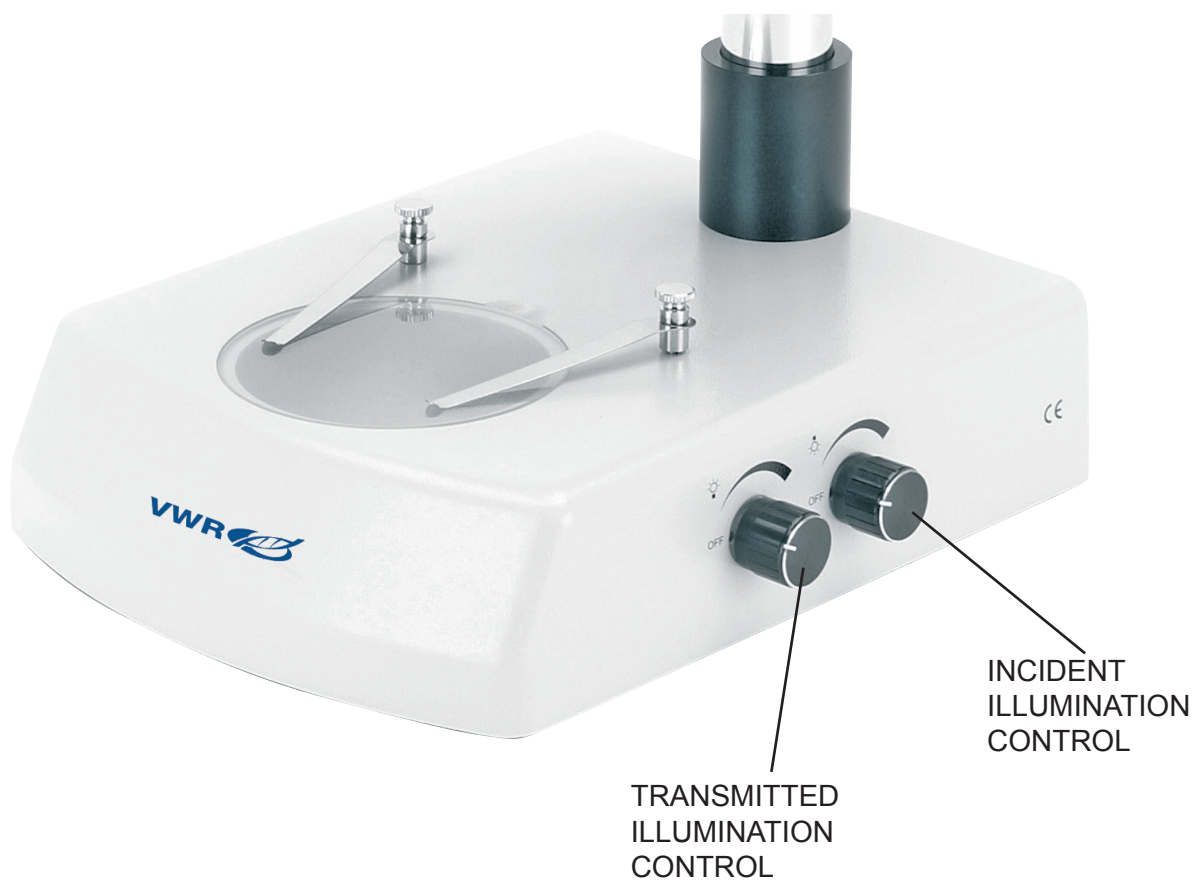Operation

Illumination system

These microscopes come with an integrated illumination system with incident and transmitted light. For the microscopes without integrated illumination systems it is possible to use an external cold light source. By using the correct light it is possible to obtain the better image of your sample.

Before starting, read the section about electrical safety precautions. Then, insert the power cord and use the brightness controls to turn on the lamps and control the light intensity.

The base is equipped with two clips that can be used to block the specimen.

Adjust interpupillary distance

Hold the right and left parts of the observation head by both hands and adjust the interpupillary distance by moving the two parts until one circle of light can be seen. If two circles appear, the interpupillary distance is too big, and if two overlapped circles appear, the interpupillary distance is too small.

Focus and regulation of the focusing tension

Put the sample to be observed on the plate and turn the zoom to the highest magnification using the bilateral zoom knob. Try to focus the sample using the focusing knobs.

Dioptric compensation

This compensation makes it possible for people with glasses to adjust the microscope to their eyes and use the microscope without glasses. Turn the zoom down to the lowest magnification. Adjust the diopter compensation ring of the right eyepiece tube until the image of the right eyepiece is clear and sharp. Repeat the procedure for the left eyepiece. Then, check the focus of the image for the whole zoom range. It should now be perfectly parfocal (focus is always maintained during the change of magnification).

Magnification and working distance

The total magnification varies from 1.75x to 180x depending on the eyepieces and auxiliary objectives used. With the bilateral zoom knob, the user can change the magnification in a factor of 6.4 with perfect parfocality and the image is always centred.

Select the desired magnification by adjusting the bilateral zoom knob. Change the eyepieces and/or add an appropriate magnifying objective lens if necessary. The total magnification used can be calculated by the following equation:

Total magnification = Eyepiece magnification x Zoom magnification x Objective lens magnification

Normal working distance for the standard configuration (1x objective lens) is 95 mm.

Video capturing (optional)

Two kinds of observation heads, binocular or trinocular, are available for the series. They can all be connected to cameras via an adaptor, for digital and analogical photo and video capturing.

Before taking a picture or filming video, pull out the light path selector lever so the light will be deflected into the photo tube. At the same time, no light will enter into the right eyepiece tube for observation. Please refer to the adaptor and camera manuals for further details.

Troubleshooting

Review the information in the table below to troubleshoot operating problems.

PROBLEM	CAUSE	SOLUTION
LIGHT DOESN'T TURN ON	Power supply not plugged.	Check that the power supply cable is well inserted on the rear of the microscope.
	Potentiometer	Rotate the brightness adjustment potentiometers and check if there is an increase in the light output.
IMAGE CANNOT BE SEEN OR IS DARK.	Optical head out of focus	Change the height of the head until you reach the correct focus.
	Brightness level is low	Rotate the brightness adjustment potentiometers.
	Photo tube inserted	If not using a camera, check that the path selection lever is fully inserted into the head (photo output disabled).
IMAGE IS UNCLEAR, BLURRED OR HAS INSUFFICIENT CONTRAST.	Objectives or filters are dirty.	Wipe them clean.
	Optical head out of focus	Change the height of the head until you reach the correct focus.
	Eyepieces set incorrectly	Check the eyepieces are set to "zero" position of dioptic compensation (rotating the ring below each eyepiece).

Repair and maintenance

Microscopy environment

This microscope is recommended to be used in a clean, dry and shock free environment with a temperature of 0-40°C and a maximum relative humidity of 85 % (non condensing). Use a dehumidifier if needed.

To think about when and after using the microscope

The microscope should always be kept vertically when moving it and be careful so that no moving parts, such as the eyepieces, fall out. Never mishandle or impose unnecessary force on the microscope.

Never attempt to service the microscope yourself.

After use, turn off the light immediately, cover the microscope with the included dust-cover, and keep it in a dry and clean place.

Electrical safety precautions

Before plugging in the power supply, make sure that the supplying voltage of your region matches with the operation voltage of the equipment and that the lamp switch is in off-position.

Users should observe all safety regulations of the region. The equipment has acquired the CE safety label. However, users do have full responsibility to use this equipment safely.

Cleaning the optics

If the optical parts need to be cleaned try first to: use compressed air. If that is not sufficient: use a soft lint-free piece of cloth with water and a mild detergent. And as a final option: use the piece of cloth moistened with a 3:7 mixture of ethanol and ether. Note: ethanol and ether are highly flammable liquids. Do not use them near a heat source, near sparks or near electric equipment. Use these chemicals in a well ventilated room.

Remember to never wipe the surface of any optical items with your hands. Fingerprints can damage the optics. Do not disassemble objectives or eyepieces in attempt to clean them.

For the best results, use the VWR cleaning kit (see catalog number below).

If you need to send the microscope to manufacturer for maintenance, please use the original packaging.

User replaceable accessories and spare parts

DESCRIPTION	QUANTITY	CAT. NO.
Eyepieces WF10x/20 mm	2	630-1779
Eyepieces WF15x/15 mm	2	630-1780
Eyepieces WF20x/10 mm	2	630-1781
Eyepiece micrometric WF10x/20 mm	1	630-1782
Additional lens 0,5x (w.d. 165 mm)	1	630-1783
Additional lens 0,75x (w.d. 117 mm)	1	630-1784
Additional lens 1,5x (w.d. 47 mm)	1	630-1786
Additional lens 2x (w.d. 26 mm)	1	630-1786
Polarising set (filters and rotating stage)	1	630-1787
Darkfield condenser	1	630-1788
Sample clip	1	630-1789
Moving stage	1	630-1790
Reflex camera adapter for full frame sensor	1	630-1791
CCD camera adapter (for 1/3" sensors)	1	630-1792
CCD camera adapter (for 1/2" sensors)	1	630-1793
Eye cups type 2	2	630-1794
White/black object-plate, type 2 dia. 95 mm	1	630-1795
Glass stage, type 2, dia. 95 mm	1	630-1796
Halogen bulb, 12V/15W	1	630-1797
Halogen bulb, 12V/15W, with dichroic mirror	1	630-1798
Dust cover type 13	1	630-1799
Heating stage, with digital temperature controller	1	630-1800
APS-C Reflex camera adapter	1	630-1801
Cleaning kit	1	630-1803

Technical service

Web Resources

Visit the VWR's website at **www.vwr.com** for:

- Complete technical service contact information
- Access to VWR's Online Catalogue, and information about accessories and related products
- Additional product information and special offers

Contact us For information or technical assistance contact your local VWR representative or visit. **www.vwr.com**.

Warranty


VWR International warrants that this product will be free from defects in material and workmanship for a period of five (5) years from date of delivery. If a defect is present, VWR will, at its option and cost, repair, replace, or refund the purchase price of this product to the customer, provided it is returned during the warranty period. This warranty does not apply if the product has been damaged by accident, abuse, misuse, or misapplication, or from ordinary wear and tear. If the required maintenance and inspection services are not performed according to the manuals and any local regulations, such warranty turns invalid, except to the extent, the defect of the product is not due to such non-performance.

Items being returned must be insured by the customer against possible damage or loss. This warranty shall be limited to the aforementioned remedies. IT IS EXPRESSLY AGREED THAT THIS WARRANTY WILL BE IN LIEU OF ALL WARRANTIES OF FITNESS AND IN LIEU OF THE WARRANTY OF MERCHANTABILITY.

Compliance with local laws and regulations

The customer is responsible for applying for and obtaining the necessary regulatory approvals or other authorizations necessary to run or use the Product in its local environment. VWR will not be held liable for any related omission or for not obtaining the required approval or authorization, unless any refusal is due to a defect of the product.

Disposal

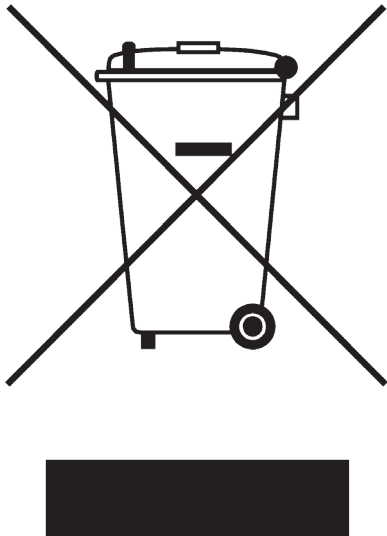This equipment is marked with the crossed out wheeled bin symbol to indicate that this equipment must not be disposed of with unsorted waste.

Instead it is your responsibility to correctly dispose of your equipment the end of its life cycle by handling it over to an authorized facility for separate collection and recycling. It is also your responsibility to decontaminate the equipment in case of biological, chemical and/or radiological contamination, so as to protect from health hazards the persons involved in the disposal and recycling of the equipment.

For more information about where you can drop off your waste equipment, please contact your local dealer from whom you originally purchased this equipment.

By doing so, you will help to conserve natural and environmental resources and you will ensure that your equipment is recycled in a manner that protects human health.

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Ireland / Northern Ireland

VWR International Ltd /
VWR International (Northern Ireland) Ltd
Orion Business Campus
Northwest Business Park
Ballycoolin
Dublin 15
Tel.: 01 88 22 222
Fax: 01 88 22 333
E-mail: sales@ie.vwr.com

Italy

VWR International PBI S.r.l.
Via San Giusto 85
20153 Milano (MI)
Tel.: 02-3320311/02-487791
Fax: 800 152999/02-40090010
E-mail: info@it.vwr.com

The Netherlands

VWR International B.V.
Postbus 8198
1005 AD Amsterdam
Tel.: 020 4808 400
Fax: 020 4808 480
E-mail: info@nl.vwr.com

New Zealand

Global Science - A VWR Company
241 Bush Road
Albany 0632, Auckland
Tel.: 0800 734 100
Fax: 0800 999 002
E-mail: sales@globalscience.co.nz

Norway

VWR International AS
Haavard Martinsens vei 30
0978 Oslo
Tel.: 02290
Fax: 815 00 940
E-mail: info@no.vwr.com

Poland

VWR International Sp. z o.o.
Limbowa 5
80-175 Gdansk
Tel.: 058 32 38 200 do 204
Fax: 058 32 38 205
E-mail: info@pl.vwr.com

Portugal

VWR International -
Material de Laboratório, Lda
Edifício Neopark
Av. Tomás Ribeiro, 43- 3 D
2790-221 Carnaxide
Tel.: 21 3600 770
Fax: 21 3600 798/9
E-mail: info@pt.vwr.com

Singapore

VWR Singapore Pte Ltd
18 Gul Drive
Singapore 629468
Tel.: +65 6505 0760
Fax: +65 6264 3780
E-mail: sales@sg.vwr.com

Spain

VWR International Eurolab S.L.
C/ Tecnología 5-17
A-7 Llinars Park
08450 - Llinars del Vallès
Barcelona
Tel.: 902 222 897
Fax: 902 430 657
E-mail: info@es.vwr.com

Sweden

VWR International AB
Fagerstagatan 18a
163 94 Stockholm
Tel.: 08 621 34 00
Fax: 08 621 34 66
E-mail: kundservice@se.vwr.com

Switzerland

VWR International GmbH
Lerzenstrasse 16/18
8953 Dietikon
Tel.: 044 745 13 13
Fax: 044 745 13 10
E-mail: info@ch.vwr.com

Turkey

Pro Lab Laboratuvar Teknolojileri Ltd.Şti.
a VWR International Company
Orta Mah. Cemal Gürsel Caddesi
Ördekcioglu İşmerkezi No.32/1
34896 Pendik - İstanbul
Tel.: +90216 598 2900
Fax: +90216 598 2907
Email: info@pro-lab.com.tr

UK

VWR International Ltd
Customer Service Centre
Hunter Boulevard - Magna Park
Lutterworth
Leicestershire
LE17 4XN
Tel.: 0800 22 33 44
Fax: 01455 55 85 86
E-mail: uksales@uk.vwr.com

**GO TO VWR.COM FOR THE
LATEST NEWS, SPECIAL OFFERS
AND DETAILS OF YOUR LOCAL
VWR DISTRIBUTOR**

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