# Cooled incubator VWR INCU-Line IL 23R Cool

## INSTRUCTION MANUAL

**European Catalogue Number:**  
**390-0881 IL 23R Cool digital**

Version: 3

Issued: 14.08.2017

CE

## Legal Address of Manufacturer

### **Europe**

VWR International bvba  
Researchpark Haasrode 2020  
Geldenaaksebaan 464  
B-3001 Leuven  
+ 32 16 385011  
<http://be.vwr.com>

Country of origin: SLOVENIA

## Table of Contents

Warning .....	4
Safety information.....	4
Package Contents .....	4
Unpacking.....	5
Installation.....	5
Intended Use .....	5
Symbols and conventions .....	5
Product specifications .....	6
Overview.....	6
Description of front panel.....	6
Instructions for use .....	7
Getting started .....	7
Operation .....	7
Temperature re-adjustment .....	7
Troubleshooting .....	8
Cleaning and maintenance .....	8
Accessories .....	8
Technical Service .....	8
Warranty .....	9
Compliance with local laws and regulations.....	9
Equipment disposal .....	9
Appendix.....	10
Your Distributor.....	12

## Warning

Configuring and operating this equipment requires ***precise knowledge*** and ***observance*** of the instructions in this ***user manual***. The equipment is designed only for its intended application.

**Important safety symbols** used in this manual:

**Warning!**

This is a warning of risk situations and dangers. Failure to observe this warning could be life-threatening. These warnings must be heeded.

This equipment is designed for 230V (50/60Hz).

Make sure it is connected to the correct power supply.

Disconnect the power cord before carrying out maintenance or cleaning work.

## Safety Information

The responsibility for operating the equipment lies exclusively with the owner or user if said equipment is incorrectly serviced, maintained or altered by persons not employed by an authorized dealer, or if the equipment is used in a manner contrary to its specified purpose.

The equipment must be maintained and operated in accordance with this user manual.

Ensure that this user manual is accessible at all times and has been read and understood.

To avoid electric shock do not spill any liquids on the equipment.

## Package Contents

Description	Quantity
Incubator IL 23R Cool	1
Stainless steel shelf	2
Power cord	1
Test certificate	1

## Unpacking

Carefully inspect the unit and accessories thoroughly upon receipt for any damage that may have occurred during shipping. Severe damage to the packaging container may indicate damage to the unit. If you suspect damage to the unit, immediately file a claim with the carrier in accordance with their instructions before contacting VWR. Also verify that all accessories are included and that the unit is in good working order before discarding the shipping packaging.

## Installation

Your satisfaction and safety require a complete understanding of this unit. Read the instructions thoroughly and be sure all operators are given adequate training before attempting to put the unit in use.

**NOTE:** This equipment must be used only for its intended application; any alterations or modifications will void your warranty and may cause injury.

## Intended use

VWR INCU-Line IL 23R Cool is a compact general purpose convection incubator, capable of uniformly cooling and heating samples from 10 °C below ambient temperature up to a maximum of 50 °C. The unit has a transparent door, so samples can be observed without opening the door, thus preventing temperature change in the chamber.

For research use only. Not intended for any animal or human therapeutic or diagnostic use.

## Symbols and conventions

The following chart is an illustrated glossary of the symbols that are used in this manual.

	<b>CAUTION</b> This symbol indicates a potential risk and alerts you to proceed with caution.
--	--

## Product Specifications

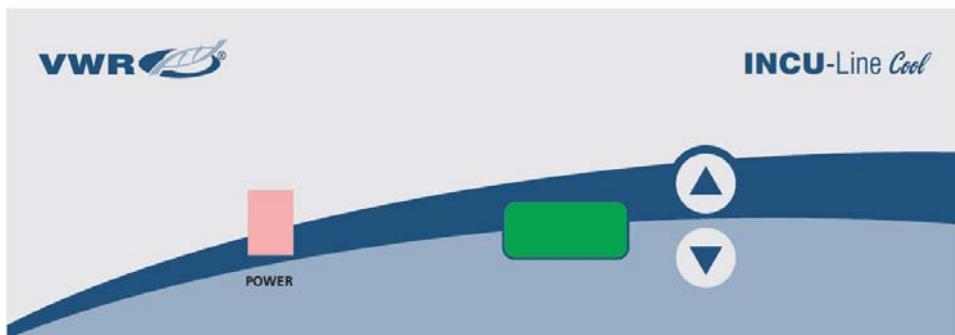
Specifications	IL 23R Cool
Electrical	230 V, 50/60 Hz, 0.4 A
Power consumption	80 W
Fuses	2 x 1 AT 250 V
Controls	PID control
Temperature range	Ambient -10 °C to +50 °C
Temperature resolution (display)	0.1 °C
Temperature accuracy at 37 °C	± 0.5 °C
Temperature fluctuation (time) at 37 °C	± 0.2 °C
Temperature variation (spatial) at 37 °C	± 1.0 °C
Environment temperature	+15 °C to +40 °C
Interior volume	23 l
Maximum load per shelf	3 kg
Maximum load per chamber	9 kg
Exterior dimensions W x D x H	380 x 470 x 450 mm
Interior dimensions W x D x H	294 x 300 x 270 mm
Weight	19.2 kg

## Overview

Access port ( $\varnothing$  7 mm) for external temperature probe

## Description of front panel

POWER - power switch: For turning the incubator ON and OFF. It lights when incubator is switched ON.

▲ - UP key: For increasing the set-temperature.

▼ - DOWN key: For decreasing the set-temperature.

Display: For displaying the actual temperature (dot blinks when you change set temperature).

## Instructions for use

### Getting Started

Check the data plate for voltage, cycle, wattage and ampere requirements. If matched to your power source, plug the power cord into a grounded outlet. Voltage should not vary more than 10 % from the data plate rating.

In selecting a location, consider all conditions which might affect performance, such as heat from radiators, autoclaves, etc. Avoid direct sun, fast-moving air currents, heating/cooling ducts and high traffic areas. Allow a minimum of 50 mm between the unit and walls or partitions which might obstruct free airflow.

### Operation

**Do not use this incubator in explosive or flammable environments. Do not heat or incubate flammable, explosive or highly reactive materials in this incubator as serious injury may result.**

Place the unit on a level surface and plug the unit into a properly grounded outlet of appropriate voltage.

Install the shelf at the desired height, with the shelf edges up. If you intend to place the samples on the bottom of the chamber, install the shelf so that it sits with the edges down on the chamber bottom, thus creating an air gap between the chamber bottom and your samples.

Place the samples in the chamber and close the door. Make sure the door is securely closed.

Turn the power switch to the ON position. The switch should illuminate. On display all "888" are displayed at first, then software version "VX.X", and then the current temperature in the chamber. Incubator starts to cool or to heat, depending on the set temperature.

Set desired temperature with UP or DOWN key. You must press and hold UP or DOWN key for more than 2 seconds, until DOT on DISPLAY starts to blink. When DOT starts to blink (it blinks for 3 seconds), set the temperature to desired value, between 10 °C below ambient temperature and up to a maximum of 50 °C. When desired temperature is set, display will automatically show the current temperature after 3 seconds (dot stops to blink).

You can always display the SET temperature by pressing and holding UP or DOWN key for more than 2 seconds.

### Temperature re-adjustment

Incubator has software for temperature re-adjustment. You must first measure the temperature in the middle of the chamber or in the sample for minimum of 2 hours. You must measure the temperature with a certificated reference temperature measuring instrument with precision 0.1 °C or more. After 2 hours or more read the temperature on thermometer and compare it with the temperature on LED display. Difference between thermometer and display is a value, which you will enter in the software.

Example 1: Reference temperature is 37.9 °C, temperature on display is 37 °C. Difference is 37.9 - 37 = 0.9  
Value is 0.9 (value which you enter in the software).

Example 2: Reference temperature is 36.2 °C, temperature on display is 37 °C. Difference is 36.2 - 37 = -0.8  
Value is - 0.8 (value which you enter in the software).

#### Procedure for temperature re-adjustment:

Press UP and DOWN key at the same time for more than 5 seconds - message "COr" (correction) appears on display and DOT starts to blink.

With UP or DOWN key set the value (when dot blinks). Pay attention on the sign of the value (plus or minus). After 3 seconds, when you set the temperature, display will automatically show the current temperature.

## Troubleshooting

Review the information in the table below to troubleshoot operating problems.

Display	Cause	Solution
Er1	Temperature probe fault, electronic board fault or incorrect setting of parameters.	Contact your distributor.
Er4	Overheating of the electronics.	Turn the main switch off and wait for the device to cool down. Then you can use the incubator again.

## Cleaning and maintenance

**Prior to any cleaning or maintenance of the unit, disconnect the power cord from the wall outlet.**

When you are using the unit for cooling, some drops of water could condense on the inside of the chamber. Use a soft cloth to wipe the inside of the chamber, if necessary.

Clean the unit internally and externally with a damp cloth and mild soap. Do not immerse the unit or spill liquids into or on the unit. Use of alcohol or ammonia-based cleaners on the door may cause it to crack. Allow the incubator to dry completely inside and outside, before reconnecting the power and using it again.

## Accessories

Description	Quantity	Cat. No.
Extra shelf, 286 x 285 mm, stainless steel	1	390-0483
Tube rack, 2 x 6 positions, Ø 32 mm	1	390-0386

## Technical service

### Web Resources

Visit the VWR's website at [www.vwr.com](http://www.vwr.com) for:

- Complete technical service contact information.
- Access to VWR's Online Catalogue, and information about accessories and related products.
- Additional product information and special offers.

### Contact us

For information or technical assistance contact your local VWR representative or visit [www.vwr.com](http://www.vwr.com).

## **Warranty**

**VWR International** warrants that this product will be free from defects in material and workmanship for a period of two (2) years from date of delivery. If a defect is present, VWR will, at its option and cost, repair, replace, or refund the purchase price of this product to the customer, provided it is returned during the warranty period. This warranty does not apply if the product has been damaged by accident, abuse, misuse, or misapplication, or from ordinary wear and tear. If the required maintenance and inspection services are not performed according to the manuals and any local regulations, such warranty turns invalid, except to the extent, the defect of the product is not due to such non-performance.

Items being returned must be insured by the customer against possible damage or loss. This warranty shall be limited to the aforementioned remedies. IT IS EXPRESSLY AGREED THAT THIS WARRANTY WILL BE IN LIEU OF ALL WARRANTIES OF FITNESS AND IN LIEU OF THE WARRANTY OF MERCHANTABILITY.

## **Compliance with local laws and regulations**

The customer is responsible for applying for and obtaining the necessary regulatory approvals or other authorizations necessary to run or use the Product in its local environment. VWR will not be held liable for any related omission or for not obtaining the required approval or authorization, unless any refusal is due to a defect of the product.

## **Equipment disposal**

This equipment is marked with the crossed out wheeled bin symbol to indicate that this equipment must not be disposed of with unsorted waste.

Instead it's your responsibility to correctly dispose of your equipment at lifecycle-end by handing it over to an authorized facility for separate collection and recycling. It's also your responsibility to decontaminate the equipment in case of biological, chemical and/or radiological contamination, so as to protect from health hazards the persons involved in the disposal and recycling of the equipment.

For more information about where you can drop off your waste of equipment, please contact your local dealer from whom you originally purchased this equipment.

By doing so, you will help to conserve natural and environmental resources and you will ensure that your equipment is recycled in a manner that protects human health.

Thank you

## Appendix

### EC Declaration of Conformity

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We, **VWR International bvba**,  
Nous, **Researchpark Haasrode 2020**,  
Wir, **Geldenaaksbaan 464**,  
**B-3001 Leuven**

hereby declare that the products, mentioned in Schedule 1 comply with the essential requirements of the following Directives of the European Parliament and council:

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and are in conformity with the following standard(s) and/or other normative document(s):

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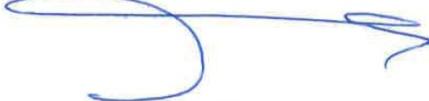
### India

VWR Lab Products Private Limited  
 No.139. BDA Industrial Suburb,  
 6th Main, Tumkur Road, Peenya Post,  
 Bangalore, India – 560058  
 Tel.: +91-80-28078400  
 Email: vwr\_india@vwr.com

### Ireland / Northern Ireland

VWR International Ltd / VWR  
 International (Northern Ireland) Ltd  
 Orion Business Campus  
 Northwest Business Park  
 Ballycoolin  
 Dublin 15  
 Tel.: +353 01 88 22 222  
 Email sales.ie@vwr.com

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VWR International GmbH
VWR International China Co., Ltd
VWR Singapore Pte Ltd